

#38 Linking m/z Protein Images with Proteomics Data

Overview

Moving forward from traditional histological and western blotting techniques, the visualization of biomolecules in tissues can now be supplemented by specific molecular identification and localization thanks to the highly sensitive discovery tool of MALDI Imaging Mass Spectrometry (MALDI IMS).

MALDI IMS consists of coating a fresh frozen or fixed tissue sample with a MALDI matrix that allows for desorption and ionization of the molecules present. A laser is then irradiated on the tissue at discrete points, thus generating an m/z spectrum across the x and y coordinates. Ion images can then be matched with an image of the same stained tissue, therefore granting valuable information of the spatial-anatomical distribution of the molecule of interest. In particular, MALDI IMS of intact proteins is of relevance, as many of the protein isoforms (proteoforms [1]) can be identified and spatially resolved.

Confirmation of protein identity remains to be one of the biggest challenges in proteomic research, with several strategies developed including bottom-up, top-down and indirect identification. The highest performance on mass resolution are provided by the Fourier transform ion cyclotron resonance (FTICR, [2]) and orbital trapping (Orbitrap, [3]) mass analyzers. However, due to their m/z range limitations, these instruments have not yet been utilized for MALDI IMS of intact proteins.

Intended Use Of This Application Note

This application note describes the utilization of MALDI FTICR MS for the generation of images of endogenous intact proteins of up to 12 KDa, with a mass accuracy of <5ppm. HTX, the manufacturers, and the users that have accepted to share their data do not guarantee the performance of the work-flow, and each lab is responsible for applying the corresponding health and safety regulations. For research purpose only.

Background

Immune cells have the capacity to attack pathogens via a wide range of strategies including reactive oxygen species (ROS) generation and chelation of metals that are essential for bacterial proliferation and antioxidant defense [4]. Calprotectin is a Zn and Mn-binding protein heterodimer, highly abundant in neutrophils, with antimicrobial and antifungal properties [5] based on the sensitization of the bacterium to superoxide oxidative damage

In the infectious foci, calprotectin is in close proximity to ROS-generating proteins, thereby being exposed to high levels of oxidative stress. Understanding the consequences of oxidative modifications on protein function of S100A8 is of high relevance to gain insights into host-pathogen interactions.

As a proof-of-concept case study, here the user was interested in imaging the various proteoforms of a particular subunit of calprotectin (S100A8), in kidney tissue from mice infected with *Staphylococcus aureus*.

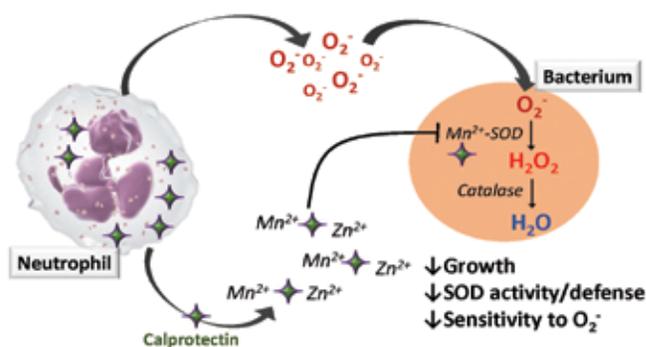


Figure 1. Adapted from [5], inhibition of bacterial processes through Mn²⁺ chelation by calprotectin. At the site of infection, neutrophils deliver a "double hit" to *S. aureus* by releasing calprotectin (crosses), which chelates Mn²⁺ and Zn²⁺, thus sensitizing *S. aureus* to ROS generated by the neutrophil.

APPLICATION NOTE

Imaging Workflow

10 μm cryosections of mouse kidney were placed onto conductive ITO coated slides (Delta Technologies). Tissue was then washed with 70% EtOH for 30 sec, 100% EtOH for 30 sec, Carnoy fluid (6:3:1 EtOH: chloroform: acetic acid) for 2 min, 100% EtOH for 30 sec, H₂O with 0.2% TFA for 30 sec, and 100% EtOH for 30 sec, and stored at -80 °C until IMS analysis.

Tissue sections were sprayed using an HTX TM-Sprayer (HTX Technologies), as described below:

| | |
|--------------------------|----------------------------|
| Flow Rate | 0.15 mL/min |
| Spray Nozzle Velocity | 1300 mm/min |
| Spray Nozzle Temperature | 30°C |
| Track Spacing | 2 mm |
| Number of Passes | 4, with offsets, rotations |
| Nitrogen Pressure | 10 psi (0.7 bar) |

Images were collected with a laser setting of $\sim 50 \mu\text{m}$ and a pixel spacing of $75 \mu\text{m}$ in x and y axis, over a mass range of m/z 1,000 to 15,000. Calibration was performed externally using CsI clusters.

Experimental Summary

| | |
|-------------------|---|
| Tissue type | Mouse kidney |
| Preservation | Mouse kidney |
| Tissue cut | 10 μm thickness |
| Matrix deposition | DHA (15 mg/mL) in 8:1:1 acetone:H ₂ O: acetic acid |
| MALDI Laser | Smartbeam II 2KHz Nd:YAG |
| Acquisition mode | Linear positive ion |

Instrumentation and Supplies

| | |
|------------------|----------------------------------|
| Cryostat | Cryostar™ Thermo NX70 |
| MALDI plate | ITO coated slides |
| Matrix | Fisher Scientific #114810250 |
| Matrix Sprayer | HTX TM-Sprayer™ |
| MALDI MS | BRUKER Solarix 15T FTICR MS |
| Imaging software | BRUKER Daltonics flexImaging 4.1 |

Results and MALDI MS Images

Fig. 2 shows MALDI FTICR IMS of intact proteins obtained from a mouse kidney infected with *S. aureus*. Data were collected with a resolving power of $\sim 75,000$ at m/z 5,000, resulting in 2552 peaks detected between m/z 2,000 and 12,000. Ions up to m/z $\sim 12,000$ were detected with high sensitivity. The average mass spectrum of the entire data set for a wild type (Wt), and a calprotectin-knock-out (KO) mouse infected with *S. aureus* are shown in Fig. 2A. The protein subunit S100A8 from the heterodimer calprotectin was detected at m/z 10,164.03 ($[M+H]^+$), -2.1 ppm) only in the Wt, and was identified by top-down fragmentation.

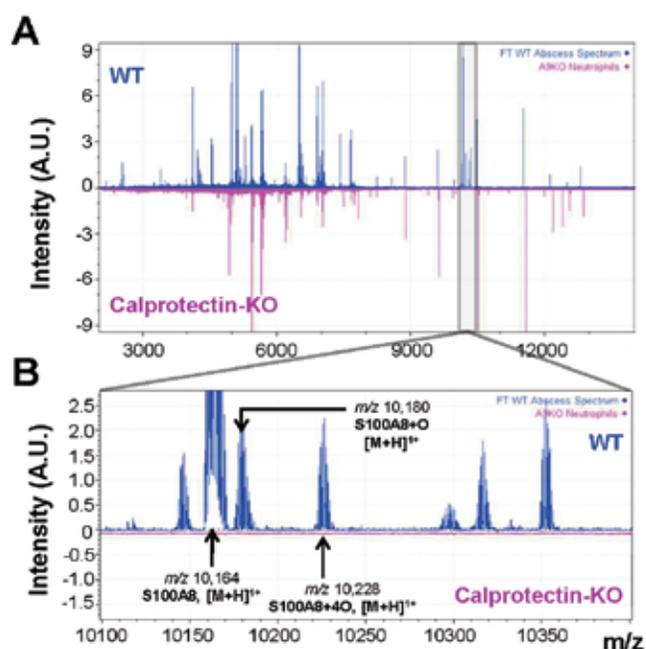


Figure 2. MALDI FTICR IMS of intact proteins from mouse kidney tissue of wild-type (blue), or calprotectin-KO (pink), infected with *S. aureus*. Images were taken from a lesion where neutrophils accumulate, in positive ion mode.

References

1. Smith, L. M., Kelleher, N. L., *The Consortium for Top Down Proteomics, Proteoform: a single term describing protein complexity.* Nat. Methods 10, 186–187 (2013)
2. Marshall, A. G., Hendrickson, C. L., Jackson, G. S.: *Fourier transform ion cyclotron resonance mass spectrometry: a primer.* Mass Spectrom. Rev. 17, 1–35 (1998)
3. Zubarev, R. A., Makarov, A.: *Orbitrap mass spectrometry.* Anal. Chem. 85, 5288–5296 (2013)
4. Russell, D.G.: *Staphylococcus and the healing power of pus.* Cell Host Microbe 3, 115–116 (2008)
5. Hood, M. I., Skaar, E. P.: *Nutritional immunity: transition metals at the pathogen-host interface.* Nat. Rev. Microbiol. 10, 525–537 (2012)
6. Gomes, L. H., Raftery, M. J., Yan, W. X., Goyette, J. D., Thomas, P.S., Geczy, C.L.: *S100A8 and S100A9-oxidant scavengers in inflammation.* Free Radic. Biol. Med. 58, 170–186 (2013)
7. Spraggins, J. M., Rizzo, D. G., Moore, J. L., Rose, K. L., Hammer, N. D., Skaar, E. P., Caprioli, R. M., *MALDI FTICR IMS of intact proteins: Using mass accuracy to link protein images with proteomics data.* J. Am. Mass Spectrom. 26, 974–985 (2015)

A closer view within the m/z range 10,160 – 10,320 allowed to identify several potential proteoforms of S100A8. Fig. 2B shows the S100A8 peak, as well as two of its oxidation products S100A8+O (m/z 10,180.07, [M+H]¹⁺), and S100A8+4O (m/z 10,228.00, [M+H]¹⁺).

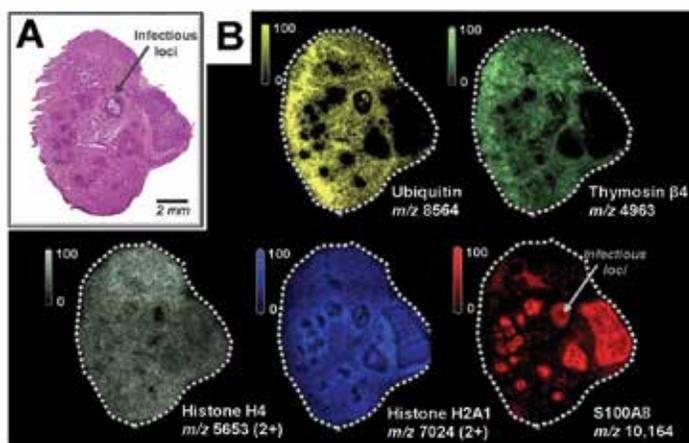


Figure 3. Selected MALDI FTICR MS ion images of intact proteins from mouse kidney infected with *S. aureus*. m/z 8,564 (ubiquitin, yellow), m/z 4,963 (thymosin 4, green), m/z 5,653 (histone H4, grey), m/z 7,024 (histone H2A1, blue), m/z 10,164 (S100A8, red). ~75 μ m spatial resolution. Acquisition time ~4 hrs. Scale bar 2 mm.

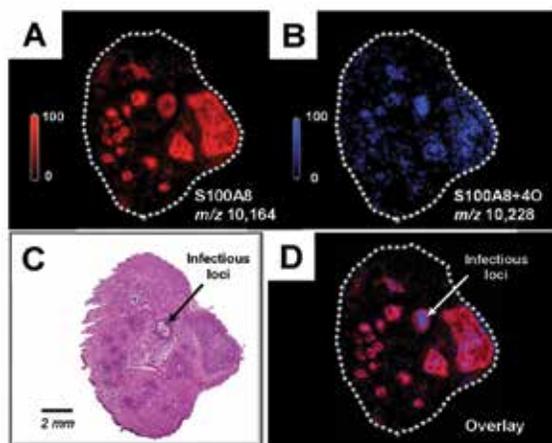


Figure 4. Selected MALDI FTICR MS ion images of (A) intact S100A8 (red, m/z 10,164) and (B) its oxidation product S100A8+4O (blue, m/z 10,228) from mouse kidney infected with *S. aureus*. (C) Shows the H&E stain, same as Fig 3A. (D) Two color overlay; m/z 10,164 (red), m/z 10,228 (blue). ~75 μ m spatial resolution. Acquisition time 2 sec / pixel, total ~4 hrs. Scale bar 2 mm.

Acknowledgements

This Technical Note, including figures and illustrations, was produced by Maria J. Torres, East Carolina Diabetes and Obesity Institute, East Carolina University, based on a consultation with Jessica Moore, co-author of the study published in [7]. Tissue images and MS data were provided courtesy of Jessica Moore, and Jeff Spraggins, Mass Spectrometry Research Center, Department of Chemistry, Vanderbilt University School of Medicine, Nashville, TN, USA.

A traditional H&E stain of the tissue is presented in Figure 3A for spatial correlation. The ion images for a series of biologically-relevant proteins (ubiquitin, thymosin, histones H4 and H2A1) and S100A8 are shown in Fig. 3B. Note the higher localization of S100A8 in the infectious loci.

Representative ion images for S100A8 and one post-translational oxidation product are shown in Fig. 4A and B, respectively. By comparing the H&E stained micrograph with the MALDI FTICR IMS overlay data (Fig. 4D), the oxidative form of S100A8 shows to accumulate near the center of the infectious foci, suggesting highly oxidative processes occur at the host-pathogen interface.

Finally, to confirm that the observed ion at m/z 10,228.00 is an oxidation proteoform of S100A8, as well as to identify the specific AA residues oxidized, top-down and bottom-up strategies were employed for identification (Fig. 5). The tryptic peptide shown in Fig 5 was bottom-up fragmented and presented oxidation of the methionine 37 and cysteine 42 residues.

MALDI FTICR IMS experiments such as the ones presented in this study are highly valuable for immunological biology given that post-translational modification of proteins (i.e. oxidation) play a key role in the immune response.


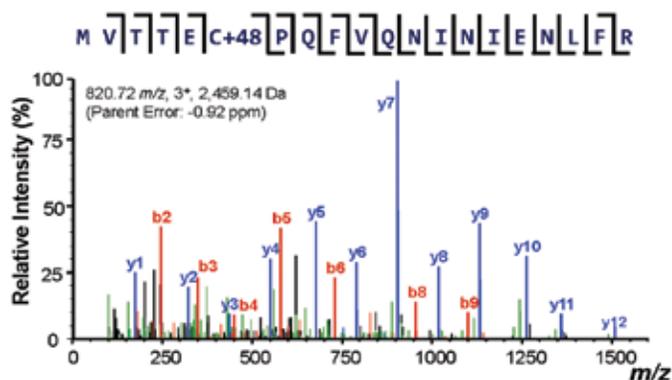


Figure 5. LC-MS/MS results for S100A8 – M370/C42O3. High-mass resolution bottom-up fragmentation data for the tryptic peptide shown above.

Conclusions

Protein identification coupled to MALDI IMS has remained a significant challenge, due to the limited mass accuracy and low resolving power. MALDI FTICR IMS significantly advances the field of molecular imaging by combining MALDI IMS and LC-MS/MS at their optimum performance.

HTX M5 Sprayer™ Tissue MALDI Sample Preparation System

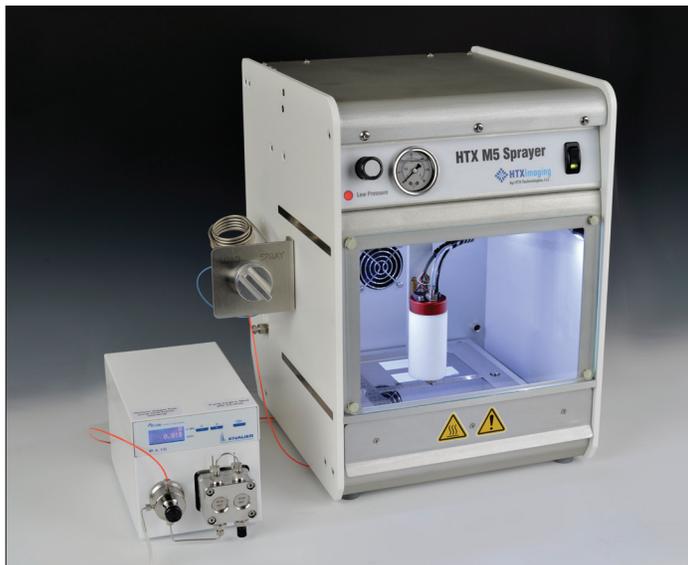

HTX M5 Sprayer™ System is an Automated MALDI Matrix Deposition System Offering **High Reproducibility** and **Superior Data Quality** for Imaging Mass Spectrometry

The HTX M5 Sprayer™ is an easy-to-use, versatile spraying system that provides automated processes for sample preparation in imaging mass spectrometry.

The proprietary spray technology of the HTX M5 Sprayer™ guarantees a very fine, uniform and consistent matrix coating crucial for high-resolution imaging and relative quantification of analytes.

The unique ability to control liquid and propulsion gas temperature creates a fine solution mist that can be deposited in a precise and adjustable pattern over all or part of any MALDI plate.

HTX M5 Sprayer™ is available worldwide exclusively from HTX Technologies, LLC.

To request further information contact:

Alain Creissen

Imaging Product Manager, HTX Technologies

acreissen@htximaging.com

HTX Technologies offers innovative sample preparation systems for advanced analytical platforms. Our integrated workflow solutions include user training, instruments, software, consumables and method development services.

Spray characteristics (wet or dry) are easily adjustable via the intuitive operator interface. Users can create and save methods for reproducible operation.

Key Characteristics

- ◆ Proprietary technology providing very small matrix droplets (<5 microns)
- ◆ High flow rate and fast sample prep (2 to 18 minutes per slide)
- ◆ Highly consistent matrix deposition across entire sample area (+/- 3% by weight)
- ◆ Unique use of temperature and nitrogen flow to control evaporation rate and matrix crystal formation
- ◆ More than 30 validated protocols covering trypsin and most matrices (e.g.: SA, CHCA, DHB, DAN, 9-AA, DHA, CMBT, THAP)
- ◆ Validated protocols for Trypsin digestion of FFPE
- ◆ Continuous matrix coverage as needed for high-resolution imaging
- ◆ Rugged operation and easy clean-up

Addressing the Matrix Deposition Challenge

The main challenge when preparing samples for MALDI Mass Spectrometry Imaging is to balance the positive effects of the matrix solution penetrating the tissue and co-crystallizing with the analyte, and the negative effects of analytes delocalization.

The all-new M5 chassis, high velocity stage and heated sample holder drawer contribute to a greater user experience and expanded process capabilities including:

- ◆ Faster and drier deposition capability
- ◆ On-tray trypsin digestion capability
- ◆ On-tray sample re-hydration



PO Box 16007 Chapel Hill, NC 27516, USA

Tel +1-919-928-5688 ◆ Fax +1-919-928-5154

info@htximaging.com ◆ www.htximaging.com